
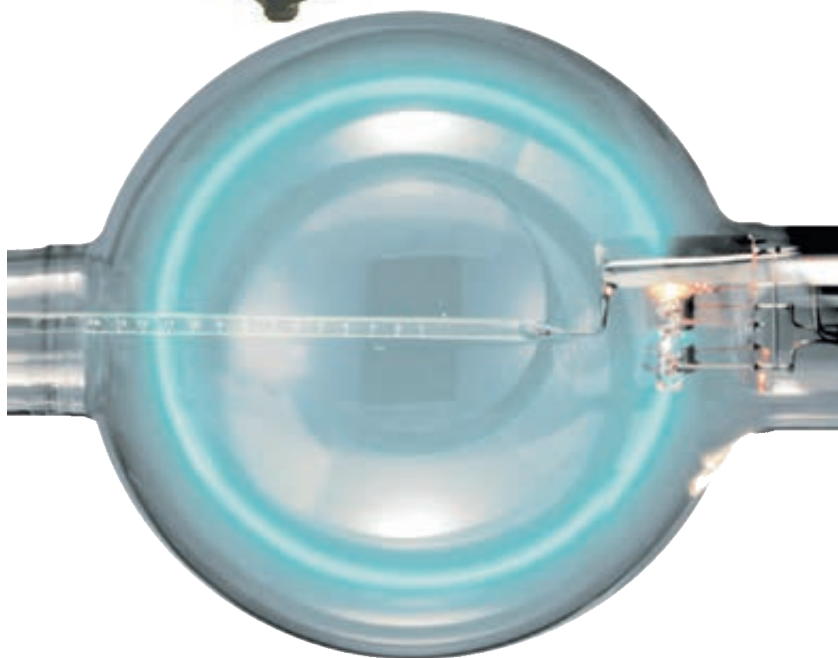


# REPLACEMENT TUBE FOR E/M APPARATUS

## INSTALLATION INSTRUCTIONS

N99-B10-7350-05

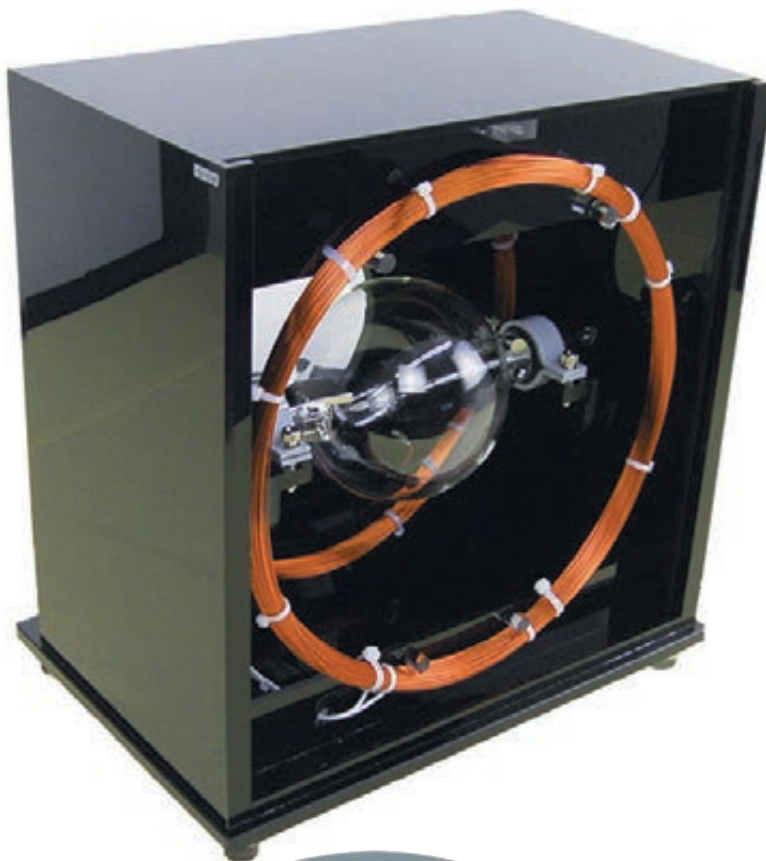

### Manual of Operations

#### **IMPORTANT!**

**Read the following before using this equipment:**

Carefully follow all instructions and observe all precautions given in this manual

Version2.1.KR072216



LOOK US UP ONLINE

# e/m TUBE INSTALLATION INSTRUCTIONS

## STEP 1


Install the discharge tube by positioning it **through the side of the Helmholtz coil** as shown in Picture 1 and 2.

**DO NOT** insert the tube through the top of the apparatus as damage may occur (as shown 3).



## STEP 2

Connect the socket to the plug of the discharge tube as shown. See the diagram to view a close up of the discharge tube connection. Pins must be positioned with holes as the socket is keyed.

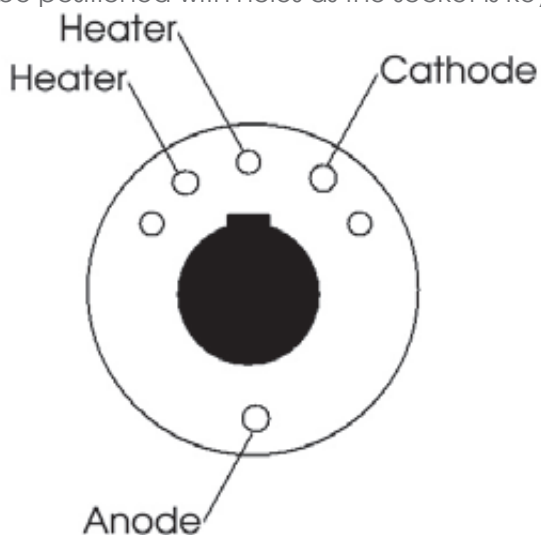

## STEP 3

Place metal fitting A (larger fitting) to the right side, and B (smaller fitting) to the left side and secure the fittings with washers and screws.

